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DE	Gebrauchshinweis – SARSTEDT Reagiergefäße	2 - 4
EN	Instructions for Use – SARSTEDT Reaction Tubes	5 - 7
BG	Инструкции за употреба – SARSTEDT реакционни епруветки	8 - 10
CS	Návod k obsluze – reagenční zkumavky SARSTEDT	11 - 13
DA	Brugsanvisning – SARSTEDT reagensbeholdere	14 - 16
EL	Οδηγίες χρήσεως – Σωληνάρια αντίδρασης SARSTEDT	17 - 19
ES	Instrucciones de uso – Microtubos SARSTEDT	20 - 22
ET	Kasutusjuhend – SARSTEDT reaktsioonikatsutid	23 - 25
FI	Käyttöohje – SARSTEDT-koeputket	26 - 28
FR	Mode d'emploi – Microtubes SARSTEDT	29 - 31
HR	Uputa za upotrebu – SARSTEDT reakcijske epruvete	32 - 34
HU	Használati utasítás – SARSTEDT reagenstartályok	35 - 37
IT	Istruzioni d'uso – Contenitori di reazione SARSTEDT	38 - 40
KO	사용 설명서 – SARSTEDT 반응 튜브	41 - 43
LT	Naudojimo instrukcijos – SARSTEDT reagentų indeliai	44 - 46
LV	Lietošanas norādes – SARSTEDT reaģentu trauki	47 - 49
NL	Gebruiksaanwijzing – SARSTEDT-reageerbuisjes	50 - 52
NO	Bruksanvisning – SARSTEDT Reagensrør	53 - 55
PL	Instrukcja obsługi – naczynia reakcyjne SARSTEDT	56 - 58
PT	Instruções de utilização – Tubos de reação SARSTEDT	59 - 61
RO	Instrucțiuni de utilizare – tuburi de reacție SARSTEDT	62 - 64
RU	Инструкция по применению – Микропробирки SARSTEDT	65 - 67
SK	Návod na Použitie – reagenčné skúmavky SARSTEDT	68 - 70
SL	Navodila za uporabo – reakcijske epruvete SARSTEDT	71 - 73
SV	Bruksanvisning – SARSTEDT mikrorör	74 - 76
TH	คำแนะนำในการใช้งาน – หลอดใส่สารทำปฏิกิริยาของ SARSTEDT	77 - 79
TR	Kullanım için talimatlar – SARSTEDT PCR Tüpleri	80 - 82
ZH	使用说明 – SARSTEDT反应管	83 - 85

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Intended use

The reaction tubes are used to hold reagents and specimen material.

They can be used for fluids such as human specimen material (e.g. nucleic acid solutions), aqueous solutions or polar solvents.

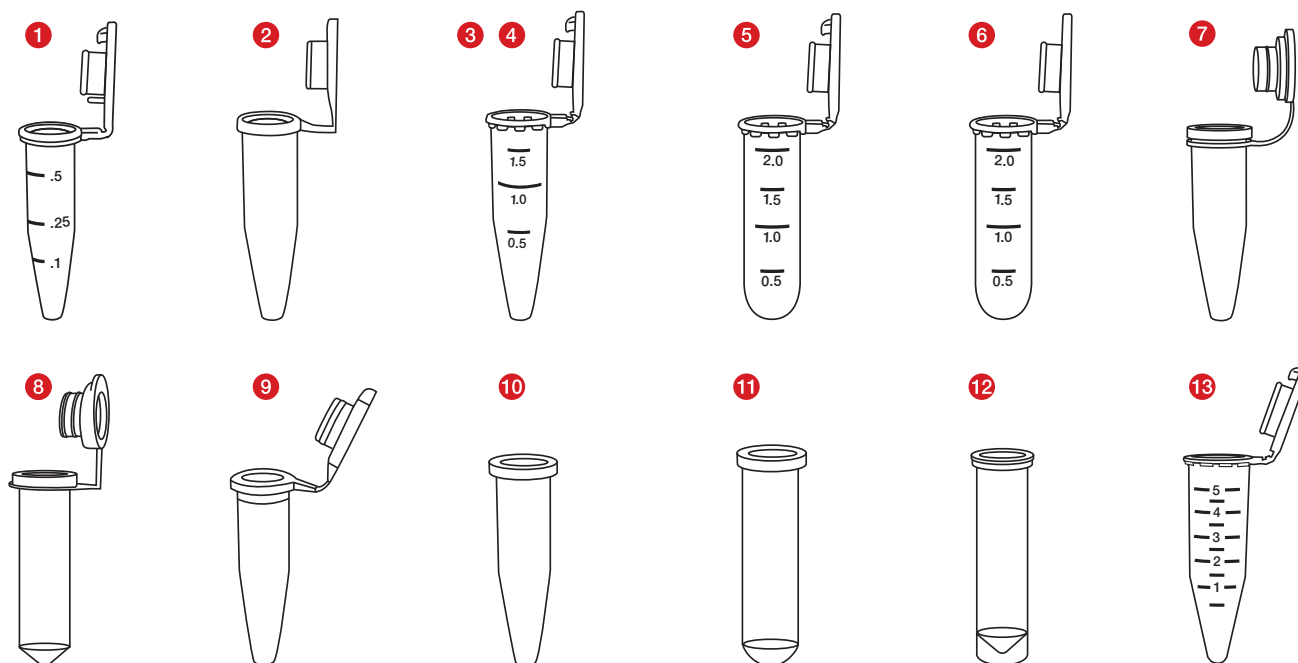
No special ambient conditions are required.

The reaction tubes should be used by qualified healthcare and laboratory personnel.

Product description


The reaction tubes are available in different variants. The designs vary in capacity and the type and properties of the cap. In addition, sterile and non-sterile variants, reaction tubes with especially low binding properties, and many different purity standards such as PCR Performance Tested and Biosphere® plus are available.

Product overview

Reaction tubes:

Type	Art. no.	Description
1	72.704.xxx	0.5 ml Safe Seal reaction tube
2	72.699.xxx	0.5 ml reaction tube
3	72.706.xxx	1.5 ml Safe Seal reaction tube
4	72.690.xxx	1.5 ml reaction tube
5	72.695.xxx	2.0 ml Safe Seal reaction tube
6	72.691.xxx	2.0 ml reaction tube
7	72.688.xxx	1.5 ml reaction tube with attached soft stopper
8	72.689.xxx	2.0 ml reaction tube with attached soft stopper
9	72.698.xxx	0.5 ml reaction tube without stopper
10	72.696.xxx	1.5 ml reaction tube without stopper
11	72.708.xxx	2.0 ml reaction tube without stopper
12	72.709.xxx	2.0 ml reaction tube, with skirted base, without stopper
13	72.701.xxx	5.0 ml reaction tube

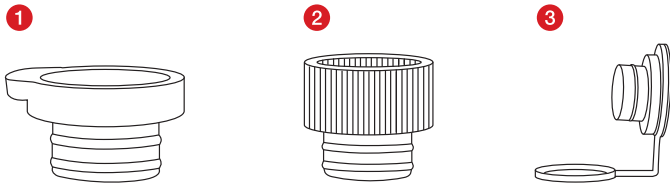Stoppers:

Type	Art. no.	Description
1	65.697.xxx	Push-in stopper, flat
2	65.723.xxx	Push-in stopper, tall
3	65.109.xxx	Soft stopper with retaining loop

Safety information and warnings

1. Personal protective equipment must be used when handling diagnostic specimens or hazardous substances.
2. Do not use reaction tubes after the expiry date.
3. Biological specimens must be handled in compliance with the applicable hygiene and safety regulations.
4. The reaction tubes are intended for single use only. After use, they must be disposed of in accordance with applicable regulations. There is a risk of contamination if reused.

Storage

Store the reaction tubes at room temperature.

Handling

1. Open the bag containing the reaction tubes and take them out.
2. Place your specimen material and any reagents into the reaction tube. Make sure that the tube is not filled above the nominal volume.
3. Close the reaction tube with the attached cap or a separate stopper.

Disposal

1. Wearing gloves and other personal protective equipment reduces the risk of infection. Follow the regulations at your facility.
2. The general hygiene guidelines and statutory regulations for the proper disposal of infectious material must be noted and complied with.
3. Contaminated reaction tubes must be disposed of in appropriate biohazard disposal containers. Follow the regulations at your facility.

Key for symbols and labels:

	Article number
	Batch number
	Use by
	CE mark
	<i>In vitro</i> diagnostic device
	Follow the instructions for use
	If reused: Risk of contamination
	Keep away from sunlight
	Store in a dry place
	Manufacturer
	Country of manufacture

Technical modifications reserved.

All serious incidents relating to the product must be reported to the manufacturer and the competent national authority.

The following is also applicable to sterile products:

	Sterilisation by irradiation
or:	
	Sterilisation by EO
	Single sterile barrier system with external protective packaging
	Do not resterilise
	Do not use if package is damaged

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