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## SARSTEDT Mikro-Schraubröhren

<b>DE</b>	Gebrauchshinweis – SARSTEDT Mikro-Schraubröhren	1 - 3
<b>EN</b>	Instructions for use – SARSTEDT Screw Cap Micro Tubes	4 - 5
<b>BG</b>	Инструкции за употреба – SARSTEDT микротръбици на винт	6 - 7
<b>CS</b>	Návod k obsluze – šroubovací mikroskopická SARSTEDT	8 - 9
<b>DA</b>	Brugsanvisning – SARSTEDT mikroreagensrør med gevind	10 - 11
<b>EL</b>	Οδηγίες χρήσεως – Μικροσωληνάρια με βιδωτό πώμα SARSTEDT	12 - 13
<b>ES</b>	Instrucciones de uso – Microtubos roscados SARSTEDT	14 - 15
<b>ET</b>	Kasutusjuhend – keermestatud mikrokatsutid SARSTEDT	16 - 17
<b>FI</b>	Käyttöohje – Kierteelliset SARSTEDT-mikroputket	18 - 19
<b>FR</b>	Mode d'emploi – SARSTEDT Microtubes à vis	20 - 21
<b>HR</b>	Uputa za upotrebu – SARSTEDT mikropruvete s navojnim grlom	22 - 23
<b>HU</b>	Használati utasítás – SARSTEDT csavaros tetejű mikrosővek	24 - 25
<b>IT</b>	Istruzioni d'uso – Microprovette con tappo a vite SARSTEDT	26 - 27
<b>KO</b>	사용 설명서 – SARSTEDT 마이크로 스크류 캡 튜브	28 - 29
<b>LT</b>	Naudojimo instrukcija – SARSTEDT mikromėgintuvėliai su užsukamais dangteliais	30 - 31
<b>LV</b>	Lietošanas norādes – SARSTEDT skrūvējamie mikrostobriņi	32 - 33
<b>NL</b>	Gebruiksaanwijzing – SARSTEDT-microbuisjes	34 - 35
<b>NO</b>	Bruksanvisning – SARSTEDT Mikrorør med skrukork	36 - 37
<b>PL</b>	Instrukcja obsługi – mikroprobówki gwintowane SARSTEDT	38 - 39
<b>PT</b>	Instruções de utilização – microtubos com tampa de rosca SARSTEDT	40 - 41
<b>RO</b>	Instrucțiuni de utilizare – micro-tuburi cu capac înșurubat SARSTEDT	42 - 43
<b>RU</b>	Инструкция по применению – Резьбовые микропробирки SARSTEDT	44 - 45
<b>SK</b>	Návod na Použitie – mikroskopická so skrutkovacím uzáverom SARSTEDT	46 - 47
<b>SL</b>	Navodila za uporabo – mikro epruveta z navojem SARSTEDT	48 - 49
<b>SV</b>	Bruksanvisning – SARSTEDT mikrorör med skruvlock	50 - 51
<b>TH</b>	คำแนะนำในการใช้งาน – หลอดเก็บตัวอย่างเกลียวขนาดเล็กของ SARSTEDT	52 - 53
<b>TR</b>	Kullanım için talimatlar – SARSTEDT Vidalı kapaklı mikro test tüpleri	54 - 55
<b>ZH</b>	使用说明 – SARSTEDT 微量螺旋管	56 - 57

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## Intended use:

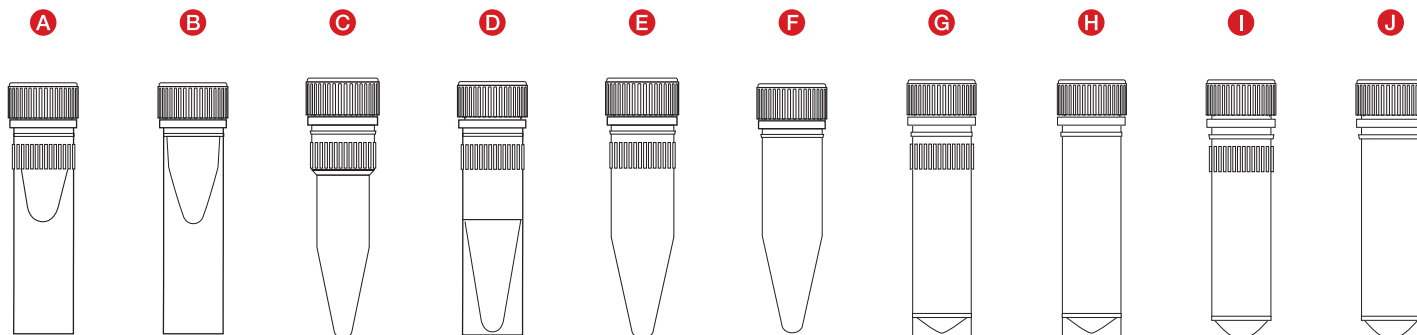
The screw cap micro tubes are used for storing reagents and specimen material. They can be used for fluids such as human specimen material (e.g. nucleic acid solutions), aqueous solutions and polar solvents.

No special ambient conditions are required.

The screw cap micro tubes should be used by qualified healthcare and laboratory personnel.


## Product description

The screw cap micro tubes are available in various designs. The different designs vary in capacity, base shape, special outer knurling and various cap options. In addition, sterile and non-sterile variants and many different purity standards such as PCR Performance Tested and Biosphere® plus are available.

## Tubes:

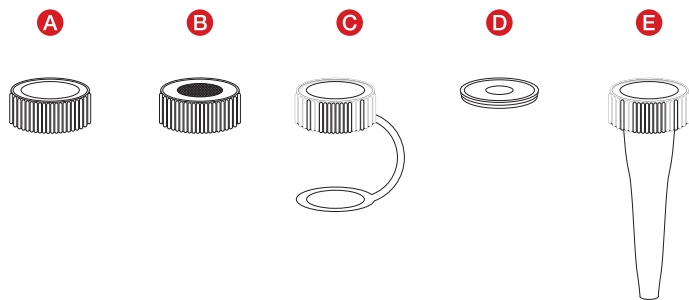Type	Art. no.	Description
A	72.730.xxx	0.5 ml screw cap micro tube, with skirted base and ridged grip
B	72.785.xxx	0.5 ml screw cap micro tube, with skirted base
C	72.733.xxx	0.5 ml screw cap micro tube, with conical base
D	72.703.xxx	1.5 ml screw cap micro tube, with skirted base and ridged grip
E	72.607.xxx / 72.692.xxx	1.5 ml screw cap micro tube, with conical base and ridged grip
F	72.687.xxx / 72.787.xxx	1.5 ml screw cap micro tube, with conical base
G	72.609.xxx / 72.694.xxx	2 ml screw cap micro tube, with skirted base and ridged grip
H	72.664.xxx	2 ml screw cap micro tube, with skirted base
I	72.608.xxx / 72.693.xxx	2 ml screw cap micro tube, with conical base and ridged grip
J	72.685.xxx	2 ml screw cap micro tube, with conical base

## Cap:

Type	Art. no.	Description
A	65.716.xxx	Screw cap for screw cap micro tubes
B	65.3716.xxx	Membrane screw cap for screw cap micro tubes
C	65.712.xxx	Attached screw cap for screw cap micro tubes
D	65.713.xxx	Colour-coded insert for attached screw cap 65.712.xxx
E	65.710.xxx	Screw cap for volume limitation

## Safety information / Warnings

1. Personal protective equipment must be used when handling diagnostic specimens or hazardous substances.
2. Do not use screw cap micro tubes after the expiry date.
3. Biological specimens must be handled in compliance with the applicable hygiene and safety regulations.
4. The screw cap micro tubes are intended for single use only. After use, they must be disposed of in accordance with applicable regulations. There is a risk of contamination if reused.

## Storage

### Storage of the screw cap micro tubes before use

Store the screw cap micro tubes at room temperature.

### Storage of the screw cap micro tubes after use

SARSTEDT screw cap micro tubes can be stored from room temperature to -80 degrees Celsius.

## Handling

1. Open the bag containing the screw cap micro tubes and take them out.
2. For variants with the cap screwed on, open the Micro screw tube by twisting the screw cap anticlockwise.
3. Place your specimen material into the screw cap micro tube. Ensure that the tube is not filled with more than the nominal volume.
4. Close the screw cap micro tube with the corresponding screw cap.
5. Store the closed screw cap micro tube according to the storage conditions specified for your specimen material.

## Disposal

1. Wearing gloves and other personal protective equipment reduces the risk of infection. Follow the regulations at your facility.
2. The general hygiene guidelines and statutory regulations for the proper disposal of infectious material must be noted and complied with.
3. Contaminated screw cap micro tubes must be disposed of in appropriate biohazard disposal containers. Follow the regulations at your facility.

## Key for symbols and labels:

	Article number
	Batch number
	Use by
	CE mark
	<i>In vitro</i> diagnostic device
	Follow the instructions for use
	If reused: Risk of contamination
	Keep away from sunlight
	Store in a dry place
	Manufacturer
	Country of manufacture

## The following is also applicable to sterile products:

	Sterilisation by irradiation
or:	
	Sterilisation by EO
	Single sterile barrier system with external protective packaging
	Sterile fluid path
	Do not resterilise

Technical modifications reserved.

All serious incidents relating to the product shall be notified to the manufacturer and the competent national authority.

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